

## Using manual syringe

1. From the DEV 2 (on rear panel) disconnect the Dual dispenser unit wire.
2. Move the XYZ-stage wire from DEV3 to DEV2
3. Remove number 1 dispenser
4. Add manual syringe holder in it's place
5. Start the tensiometer
6. Start the OneAttension software
7. Before using the syringe, do a new CALIBRATION

when done with manual syringe:

8. Close the software
9. Turn off the tensiometer
10. Remove manual syringe holder
11. Add number 1 dispenser back to it's place
12. Move the XYZ-stage wire from DEV2 to DEV3
13. Reconnect the Dual dispenser unit wire to DEV2
14. Start the tensiometer
15. Start the OneAttension software
16. Do a new CALIBRATION